



## LOG COLLECTOR HOW TO


Copy the file InstallerLogToolv114ENSEGv3.zip to C:\LabSolutions

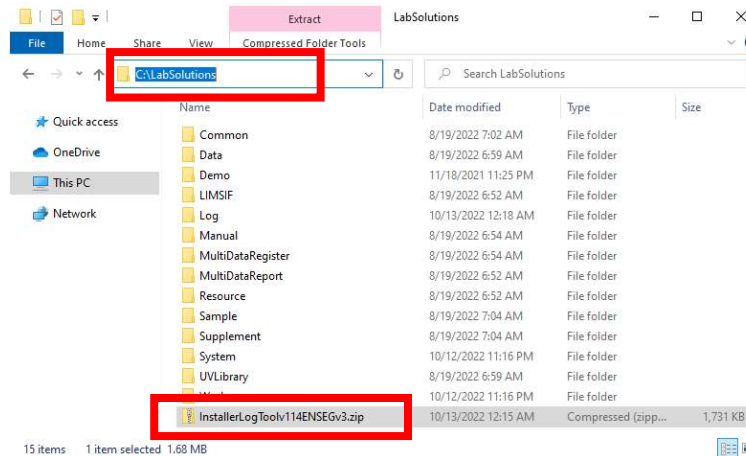

Right click on the file “InstallerLogTool114ENGSEGV3.zip” and click “Extract All...”

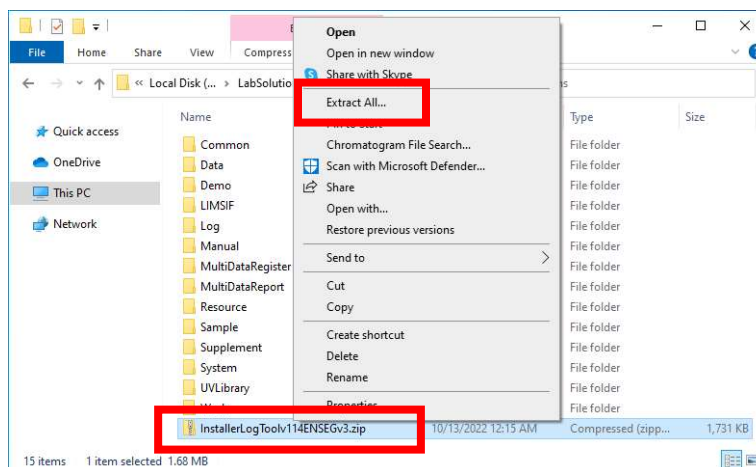


Click “Extract”

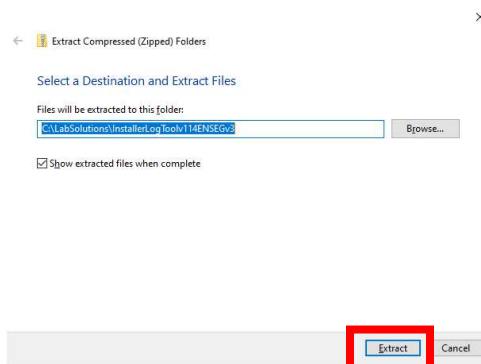


Right click on the file InstallerLogCollectorEN.bat and click “Run as administrator”

Note: If you run the tool without administrator privileges, it will terminate. In this case, no logs are collected!

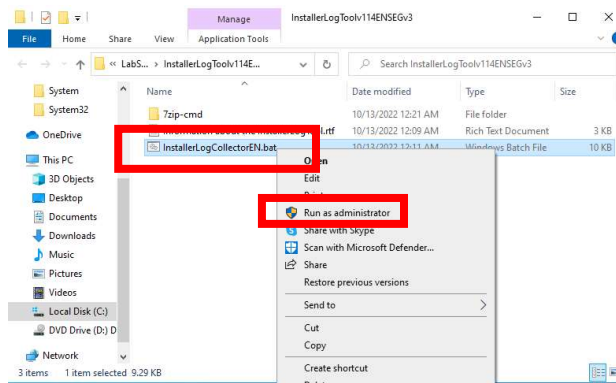

When successful a Log\_date-time-PCName.zip file is generated.

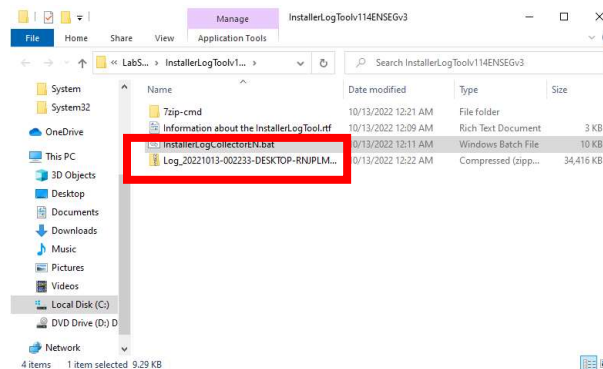

Send the Log\_date-time-PCName.zip to SEG via the provided upload link. If no upload link is provided, ask for one.